



MARS 6

Microwave Acid Digestion

Method Note Compendium

October 1, 2019

Contents

Agriculture

Alfalfa	11
Boric Acid HF Neutralization	12
Cannabis (Plant)	13
Coffee Beans	14
Cotton	15
Feed Grain	16
Fertilizer - AOAC Method 2017.02	17
Fertilizer -AOAC Method 2006.03	18
Geranium	19
Impatiens	20
Leather	21
Palm Leaves	22
Plant Tissue	23
Soybean	24
Soybean Meal	25
Spinach Leaves	26
Straw	27
Tea Extract	28
Tea Tannin	29
Tobacco	30
Wheat	31
Wood	32
Wood Pulp	33
Wool	34
Yeast	35

Clinical & Biological

Animal Tissue (Dry)	37
Animal Tissue (Wet)	38
Beef Bone	39

Blood (Human)	40
Bovine Blood	42
Bovine Liver (Wet)	43
Brain Tissue (Dog)	44
Dog Feces (Dry)	45
Fingernails	46
Hair	47
Lobster Hepatopancreas (Tort-CRM)	48
Rat Kidney	49
Urine - Human	50

Consumer Products

Boric Acid HF Neutralization	52
Candle Fragrance in iPrep	53
Cosmetics (Liquid Make-Up)	54
Cotton	55
Deodorant (Powder)	56
Eye Shadow	57
Foam (Mattress) in iPrep	58
Sanitizers (Hand)	59
Shampoo (Dandruff)	60
Soaps (Lotion and Foam)	61
Sunscreen in iPrep	62
Suppository (Capsule)	63
Toothpaste	64
Vaseline	65
Wax (Candle)	66

Environmental & Regulatory

Boric Acid HF Neutralization	68
Fertilizer - AOAC Method 2017.02	69
Fertilizer - AOAC Method 2006.03	70
Fly Ash in iPrep	71
Limestone	73

NPDES (Waste Water)	74
Pine Needles	75
RoHS (For Pb, Hg, and Cd Analysis)	76
Sediment (BCCS - 1 CRM) (Leach)	77
Sediment (Buffalo River) (Leach)	78
Sludge (Industrial)	79
Soil (Montana - CRM) (Leach)	80
TCLP Extract	81
US EPA 3015 (Aqueous Samples)	82
US EPA 3015a (Aqueous Sample)	83
US EPA 3051 (Solid Sample)	84
US EPA 3051a (Solid Sample)	85
US EPA 3052 (Sludge, Soil, Sediment or Oil)	86
USP 232/233 (Pharmaceuticals)	87
Waste Activated Sludge	88
Water (For Analysis of Phosphorous)	89
Wood	90

Cherry	108
Chewing Gum	109
Chicken (Boneless)	110
Chili Candy (Mexican)	111
Chili Powder in iPrep	112
Chips (Potato)	113
Citrus Leaves	114
Clam Puree	115
Cocoa	116
Coconut	117
Coffee Beans	118
Cucumber	119
Dog Food	120
Fish Tissue	121
Flounder	122
Flour	123
Food	124
Food Coloring (Liquid)	125
Fruit Juice	126
Gelatin	127
Granola Bar	128
Grape	129
Grapefruit	130
Gummy Bear	131
Ham	132
Infant Cereal (Rice Grain)	133
Infant Formula (Liquid) in iPrep	134
Infant Formula (Liquid)	135
Infant Formula (Powder) in iPrep	136
Infant Formula (Powder)	137
Kidney Bean	138
Kielbasa Sausage	139
Maple Syrup	140
Mayonnaise	141
Meal Replacement Shake (Chocolate)	142
Milk (Whole - Liquid)	143

Food Testing

Alum	92
Apple	93
Asparagus	94
Aspartame in iPrep	95
Bacon (Cooked)	96
Beef (Ground)	97
Beef Bouillon	98
Beer (Light)	99
Blueberry	100
Broccoli	101
Caramel Color	102
Carrot	103
Celery	104
Cheese (Cracker)	105
Cheese (Powder)	106
Cheese (Processed)	107

Milk (Powder)	144
Nutritional Drink (Adult)	145
Olive Oil	146
Orange Juice	147
Oyster Puree	148
Peanut Butter	149
Pear	150
Pepper (Bell, Chili, Etc.)	151
Pistachio Nuts	152
Pizza (Frozen, Pepperoni)	153
Plum	154
Pork (Ground)	155
Potato	156
Potato Chips	157
Pretzel (Salted)	158
Protein Bar	159
Raspberry	160
Rice	161
Safflower Oil	162
Salad Dressing (Ranch)	163
Salami	164
Sausage	165
Soda (Diet)	166
Sodium Free - Seasoning	167
Soybean	168
Spinach	169
Spinach Leaves	170
Spinach Pasta	171
Strawberry	172
Sugar (Granulated)	173
Tea (Chai) in iPrep	174
Tea Leaves	175
Tomato Leaves	176
Tomato Paste (70% Moisture)	177
Tomato Soup	178
Walnut	179

Watermelon	180
Wheat Crackers	181
Whey	182
Whey (Powder)	183
Wine	184
Yogurt (Plain)	185

Geoscience & Mining

Alpha - Alumina in iPrep	187
Alumina Beads (with Silver Substrate) in iPrep	188
Alumina Powder in iPrep	190
Aluminum Nitride	191
Aluminum Oxide	192
Aluminum Oxide (H_3PO_4) in iPrep	193
Asbestos	194
Automotive Catalyst	195
Barium Titanate ($BaTiO_3$)	196
Bastnaesite Ore	197
Bauxite	199
Bismuth Aluminate Mixture	201
Bismuth Yttrium Iron Oxide in iPrep	202
Boiler Scale (80-90% Fe)	204
Boric Acid HF Neutralization	205
Boron Carbide (B_4C) in iPrep	206
Carbon	207
Cement	208
Ceramic (Fused Silica)	209
Ceramics	210
Char - Sulfuric Acid	211
Chromium Carbide	212
Chromium Oxide (Cr_2O_3) in iPrep	213
Clay	214
Coal in iPrep	215
Coal	216
Coal Ash	217

Cryolite / Electrolytic Bath	218
Feldspar in iPrep	220
Ferric Oxide in iPrep	222
Furnace Slag	223
Gadolinium Zirconium Oxide in iPrep	224
Gas Atomized Powder in iPrep	226
Glass Fiber	227
Hafnium Oxide	228
Ilmenite (FeTiO ₃)	229
Iron Ore	230
Kaolin	231
Kerosene	232
Lead Sulfide (PbS)	233
Lithium Nickel Cobalt Manganese Oxide	234
Magnesium Oxide (MgO 40-54%)	235
Manganous Oxide (MnO)	236
Mica in iPrep	238
Mill Tailings	239
Mineral Ore	241
Molybdenum Disulfide (MoS ₂)	243
Nickel (II) Hydroxide	245
Niobium Pentoxide	246
Petroleum Coke	247
Phosphate (21%)	249
Phosphate Rock	250
Pollucite Ore in iPrep	251
Pyrophyllite (AlSi ₂) ₅ (OH)	253
Quartz Powder	254
Rock (Pulverized)	255
Rock, High Sulfide in iPrep	256
Rutile Ore	258
Sand	259
Shale in iPrep	260
Silica Sand	261
Silica Sand in iPrep	262
Silicon Dioxide	263

Silicon Nitride in iPrep	264
Silicon Wafer	265
Slag - Furnace	266
Sodium Antimonate in iPrep	267
Spodumene (Lithium Feldspar)	268
Sulfide Bulk Concentrate in iPrep	270
Tantalum Pentoxide (Ta ₂ O ₅) in iPrep	272
Titanium Dioxide	273
Titanium Slag Ore in iPrep	274
Tungsten Carbide	275
Tungsten Ore (50% W; 10% Fe; 2% Mn)	276
Tungsten Oxide	277
Vermiculite	278
Wollastonite (Calcium Silicate) in iPrep	279
Zeolite	280
Zinc Oxide (ZnO)	281
Zinc Oxide (ZnO) in iPrep	282
Zircon in iPrep	283

Industrial Hygiene

Asbestos	285
Filter (MCE)	286
Filter Paper - 37 mm Mixed Cellulose Ester	287
Filter Paper - 47 mm Polycarbonate	288
Filter Paper - Cellulose	289
Ghost Wipe	290
Paint Chips (For Pb Analysis)	291
Quartz Fiber Filters	292
Toray Filter in iPrep	293
Toray Filter (Anion)	294
Toray Filter (Cation)	295

Materials Science

Acrylic Tape in iPrep	297
-----------------------	-----

Bismuth Aluminate Mixture	298
Bismuth Yttrium Iron Oxide in iPrep	299
Ceramic Powder	301
Ceramic Powder in iPrep	302
Circuit Board (Cryo - Ground)	303
Ferric Oxide in iPrep	304
Gadolinium Zirconium Oxide in iPrep	305
Gas Atomized Powder in iPrep	307
Glass (Ground)	308
Graphite Composite - Cyanate Ester Resin	309
Graphite Composite - Epoxy Resin	310
Graphite Composite - PEEK Resin	311
Graphite Fiber - Cyanate Siloxane Resin	312
Graphite Fiber - Epoxy Resin (Fiber Content)	313
Pollucite Ore in iPrep	314
Siloxanes	316

Metals & Alloys

Alpha - Alumina in iPrep	318
Aluminum (Metal)	319
Aluminum Alloy	320
Aluminum Alloy (20% SiC)	321
Aluminum Alloy (BCS 182)	322
Aluminum Oxide	323
Aluminum Oxide (H ₃ PO ₄) in iPrep	324
Boron	325
Boron Carbide (B ₄ C) in iPrep	326
Carbonyl Iron in iPrep	327
Cobalt - Chromium Alloy (Co-Cr)	328
Copper (Metal)	330
Copper Ore	331
Gallium	332
Iron - Chromium Alloy	333
Iron - Chromium Alloy in iPrep	335
Iron - Lignosulfonate	336

Iron - Manganese Alloy	337
Iron - Silicon Alloy	338
Lead Sulfide (PbS)	339
Lithium Nickel Cobalt Manganese Oxide	340
Magnesium Oxide (MgO 40-54%)	341
Manganous Oxide (MnO)	342
Metal Alloy - FeCr	344
Metal Alloy - NiCr	345
Metal Alloy (Ag, Au, Cu) in iPrep	346
Metal Alloy (Ag, Co, Cu) in iPrep	347
Metal Alloy (Ag, Cu, Sn) in iPrep	348
Metal Ore - Precious	349
Metals Alloy (Co-Cr)	351
Nickel (II) Hydroxide	353
Nickel (Metal)	354
Niobium - Titanium Alloy	355
Phosphate (21%)	356
Platinum Metal	357
Platinum Ruthenium in iPrep	358
Rhodium Sponge in iPrep	359
Selenium Alloy	360
Sn - Ag Solder	361
Stainless and Low Carbon Steel in iPrep	362
Stainless Steel	363
Steel (Stainless)	364
Strontium Ferrite in iPrep	365
Titanium Alloy	366
Titanium Dioxide	367
Tungsten Ore (50% W; 10% Fe; 2% Mn)	368
Zinc Oxide (ZnO)	369
Zinc Oxide (ZnO) in iPrep	370
Zircon Alloy	371

Nutraceuticals

Biotin	373
--------	-----

Cannabis (Plant)	374
Chili Powder in iPrep	375
Citrus Leaves	376
Ginkgo (Ground)	377
Ginkgo (Whole Pill)	378
Infant Formula (Liquid) in iPrep	379
Infant Formula (Liquid)	380
Infant Formula (Powder) in iPrep	381
Infant Formula (Powder)	382
Kelp	383
Milk (Whole - Liquid)	384
Milk (Powder)	385
Plant Tissue	386
USP 232/233 (Pharmaceuticals)	387
Wheat	388
Whey (Powder)	389

Organic Chemicals

Di-(2-ethylhexyl)phosphoric Acid in iPrep	391
Diazo Sulfonyl Chloride	392
Food Coloring (Erythrosine Based)	393
Methylcellulose (Hydroxypropyl)	394
Monoethanolamine	395
Photoresist	396
Polyglutaramide	397

Paints & Coatings

Paint (Latex Based Liquid)	399
Paint Chips (For Pb Analysis)	400
Photoresist	401
Polyurethane	402
Polyurethane in iPrep	403
Toner (Ink) in iPrep	404

Toner (Powder) in iPrep	405
-------------------------	-----

Paper & Pulp

Filter Paper - 37 mm Mixed Cellulose Ester	407
Filter Paper - 47 mm Polycarbonate	408
Filter Paper - Cellulose	409
Paper Pulp	410

Personal Care

Boric Acid HF Neutralization	412
Cosmetics (Liquid Make-Up)	413
Eye Shadow	414
Shampoo (Dandruff)	415

Pharmaceutical & Biotech

Allergy Pill (Ground)	417
Allergy Pill (Whole Pill)	418
Antioxidant (Ground)	419
Antioxidant (Whole Pill)	420
APIs with Aromatic Ring Structures in iPrep	421
Ascorbic Acid	422
Aspirin (Ground)	423
Aspirin (Whole)	424
Beta Carotene	425
Biotin	426
Calcium Carbonate	427
Calpan Pantothenic Acid	428
Cephalexin	429
Chromium Chelate	430
Cold & Flu Medicine (Liquid)	431
Cupric Sulfate	432
Diclofenac K	433
Empty Capsule (Gel)	434

Erythromycin	435
Fish Oil (No Capsule)	436
Flurbiprofen	437
Folic Acid	438
Ginkgo (Ground)	439
Ginkgo (Whole Pill)	440
Hydrocodone	441
Ibuprofen (Ground)	442
Kelp	443
Losartan K	444
Magnesium Carbonate	445
Magnesium Oxide	446
Manganese Carbonate	447
Medroxyprogesterone	448
Metformin (Ground)	449
Mineral Oil	450
Multivitamin (Ground)	451
Multivitamin (Whole)	452
Naproxen Sodium	453
Niacinamide	454
Pantothenic Acid	455
Potassium Chelate	456
Protopic Ointment	457
Pyridoxine	458
Riboflavin	459
Selenium Chelate	460
Sleep Aid	461
Sodium Polystyrene Sulfonate in iPrep	462
Stearic Acid	463
Sudan Orange (4-Fluorophenyl) in iPrep	464
Thiamine	465
Transdermal Patch	466
Trixie Phosphate in iPrep	467
USP 232/233 (Pharmaceuticals)	468

Vitamin B-12	469
Vitamin C (Ground)	470
Vitamin C (Whole)	471
Vitamin Concentrate (High Fe)	472
Vitamin D	473
Vitamin D3	474
Vitamin E	475
Vitamins	476
Zinc Sulfate	477

Plastics, Polymers, & Oils

Acrylamide	479
Brominated Flame Retardant in iPrep	480
Bunker Oil in iPrep	481
Char - Sulfuric Acid	482
Crude Oil	483
Diesel Fuel	484
Epoxy Hardener	485
Fatty Alcohol	486
Fluorinated Ethylene Propylene (FEP) in iPrep	487
Fuel Oil (Slurry) in iPrep	488
Gasoline	490
Glycerol Ester of Wood Rosin (Ground) in iPrep	491
HDPE	492
Kerosene	493
Kevlar in iPrep	494
Lube Oil	495
Methyl Ester of Hydrogenated Wood Rosin (Ground) in iPrep	496
Mineral Oil	497
Motor Oil (New)	498
Motor Oil (Waste)	499
Motor Oil (Waste) in iPrep	500
Nylon	501
Oil Wear Metals	502
PET (Polyethylene Terephthalate)	503

PET (Polyethylene Terephthalate) in iPrep	504
Petrolatum	505
Plasticizer (Filter) in iPrep	506
Plasticizers in iPrep	507
Polycarbonate Resin	508
Polyethersulfone (Filter Paper)	510
Polyethylene - High Density	511
Polyphenylene Sulfide	512
Polypropylene	513
Polyterpene Resin in iPrep	514
Polyurethane	515
Polyurethane in iPrep	516
PVC (Polyvinyl Chloride Resin)	517
Rubber	518
Rubber (Synthetic)	519
Silane Terminated Polyether in iPrep	520
Silicone Tubing - TiO ₂ Filler	521
Styrene Acrylate	522
Terephthalic Acid	523
Wax (Petroleum Based)	524

Water & Wastewater Treatment

NPDES (Waste Water)	526
Sludge (Industrial)	527
TCLP Extract	528
US EPA 3015 (Aqueous Samples)	529
US EPA 3015a (Aqueous Sample)	530
Waste Activated Sludge	531
Water (For Analysis of Phosphorous)	532