



Discover[®] SP-D 80

Automated Microwave Digestion System

Method Note Compendium

Updated May 7, 2019

Contents

Agriculture

Alfalfa	2
Boric Acid HF Neutralization	3
Tobacco	4
Wheat	5
Whey	6
Yeast	7

Clinical & Biological

Boric Acid HF Neutralization	9
Bovine Liver	10
Lobster Hepatopancreas (Tort-2 CRM)	11

Consumer Products

Boric Acid HF Neutralization	13
Cotton	14
Deodorant (Powder)	15
Eye Shadow	16
Shampoo	17
Suppository	18
Toothpaste	19
Vaseline	20
Wax (Candle)	21

Environmental & Regulatory

Boric Acid HF Neutralization	23
Fertilizer - AOAC 2006.03	24
NPDES	25
Phosphorous	26

Pine Needles	27
ROHS	28
Sediment - Buffalo River (Leach)	29
Sediment (Leach)	30
Sludge	31
Soil - Montana (Leach)	32
TCLP	33
US EPA 3015a (Aqueous Sample)	34
US EPA 3051 (Solid Sample)	35
US EPA 3052 (Sludge, Soil, Sediment, or Oil)	36
Wood	37

Food Testing

Apple	39
Asparagus	40
Bacon	41
Beer (Light)	42
Blueberry	43
Boost Drink	44
Boric Acid HF Neutralization	45
Broccoli	46
Caramel Color	47
Carrot	48
Celery	49
Cereal (Krispies)	50
Cheese	51
Cherry	52
Chicken (Boneless)	53
Chili Candy (Mexican)	54
Chips (Potato)	55
Coffee	56
Corn Meal	57

Cracker (Saltine)	58
Cucumber	59
Dog Food (Hard)	60
Dressing (Ranch)	61
Flour	62
Food Coloring	63
Fruit Juice	64
Grape	65
Grapefruit	66
Ground Beef	67
Ham	68
Infant Formula (Powder)	69
Kidney Bean	70
Mayonnaise	71
Milk - Liquid	72
Milk - Powder	73
Olive Oil	74
Orange	75
Orange Juice	76
Peanut Butter	77
Pear	78
Pepper (Green, Bell)	79
Pizza (Frozen, Ground)	80
Plum	81
Pork (Raw)	82
Pork Gelatin	83
Potato	84
Pretzel (Salted)	85
Raspberry	86
Rice (White)	87
Safflower Oil	88
Soda (Diet)	89
Soybean	90
Spinach	91
Strawberry	92
Sugar	93

Tea Leaves	94
Tomato Leaves	95
Tomato Soup	96
Walnut	97
Watermelon	98
Wine	99
Yeast	100
Yogurt (Plain)	101

Geoscience & Mining

Boric Acid HF Neutralization	103
Coal	104
Rock	105

Industrial Hygiene

Boric Acid HF Neutralization	107
Filter Paper - 37 mm Mixed Cellulose Ester	108
Filter Paper - Polyethersulfone	109
Filter - 47 mm Polycarbonate	110
Ghost Wipe	111
Glass	112

Materials Science

Boric Acid HF Neutralization	114
Carbon Nanotubes	115
Circuit Board	116
Glass	117

Metals & Alloys

Aluminium	119
Aluminium Oxide	120
Boric Acid HF Neutralization	121

Copper	122
Nickel	123
Steel (Stainless)	124
Titanium Dioxide	125
Zinc Oxide	126

Nutraceuticals

Banana Leaves	128
Boric Acid HF Neutralization	129
Citrus Leaves	130
Corn Meal	131
Ginkgo (Ground)	132
Ginkgo (Whole)	133
Palm Leaves	134
Pine Needles	135
Pine Straw	136
Rice Flour	137
Tea Leaves	138
Tomato Leaves	139

Organic Chemicals

Boric Acid HF Neutralization	141
Photoresist	142

Paints & Coatings

Boric Acid HF Neutralization	144
Photoresist	145

Paper & Pulp

Boric Acid HF Neutralization	147
Filter Paper - 37 mm Mixed Cellulose Ester	148
Filter Paper - Polyethersulfone	149
Filter - 47 mm Polycarbonate	150
Ghost Wipe	151

Pulp (Paper)	152
--------------	-----

Pharmaceutical & Biotech

Allergy Pill (Ground)	154
Allergy Pill (Whole)	155
Antioxidant (Ground)	156
Antioxidant (Whole)	157
Aspirin (Ground)	158
Aspirin (Whole)	159
Boric Acid HF Neutralization	160
Calpan Pantothenic Acid	161
Capsule (Empty)	162
Cephalexin	163
Chromium Chelate	164
Cold and Flu Medicine (Liquid)	165
Diclofenac K	166
Fish Oil (No Capsule)	167
Folic Acid	168
Ginko (Ground)	169
Ginko (Whole)	170
Ibuprofen (Ground)	171
Kelp	172
Losataran K	173
Magnesium Oxide	174
Manganese Carbonate	175
Metformin	176
Multivitamin (Ground)	177
Multivitamin (Whole)	178
Stearic Acid	179
Thiamine	180
USP 232/233	181
Vitamin C (Ground)	182
Vitamin C (Whole)	183
Vitamin D3 (Ground)	184
Zinc Sulfate	185

Plastics, Polymers, & Oils

Acrylamide	187
Boric Acid HF Neutralization	188
Bunker Oil	189
Char - Sulfuric Acid	190
Crude Oil	191
Epoxy Hardener	192
Fatty Alcohol	193
Gasoline	194
HPDE	195
Lube Oil	196
Mineral Oil	197
Motor Oil (New)	198
Nylon	199
PET	200
Plastic	201
Polycarbonate Resin	202
Polypropylene	203
Polyurethane	204
Rubber (Synthetic)	205
Rubber (Tire)	206
Terephthalic Acid	207

Water & Wastewater

NPDES	209
TCLP	210